

# True UHV — Low Power

The CS202\*B is a True UHV cold head (10<sup>-11</sup> Torr) where all of the rubber o-ring seals have been replaced with welded joints and metal seals. A CF flange is directly welded to the cryocooler.

These True UHV systems are bakeable to 80C, and the 10K CS202AB is bakeable to 1200C if the displacer is removed. (Special Training Required).

Cold tip extensions are available to put the sample right where it is needed in the chamber. The extra cooling power of the CS204\*B comes in handy when working with long cold tip extensions.

#### **Applications**

- UHV
- Surface Science
- UHV Manipulator for XYZ motion
- Photoemission Spectroscopy

#### **Features**

- True UHV (10-11 Torr)
- Bakeable to 80C (10K version can remove displacer and bake to 200C)
- Open Sample Space
- Optional Cold Tip Extensions
- Cryogen Free
- Operation in Any Orientation
- Fully Customizable

#### **Typical Configuration**

- Cold head (DE-202AB)
- Compressor (ARS-2HW)
- 2 Helium Hoses
- Tue UHV welded stainless steel instrumentation skirt with 6" rotatable CF flange
- Nickel Plated OFHC radiation shield terminating 0.125" short of the cold tip
- 2 High purity quartz windows
- Instrumentation for temperature measurement and control:

10 pin UHV feed through

36 ohm thermofoil heater

Silicon diode sensor curve matched to (±0.5K) for control

Calibrated silicon diode sensor ( $\pm 12$  mk) with 4 in. free length for accurate sample measurement.

- Wiring for electrical experiments:
  - 10 pin hermetic feed through
  - 4 copper wires
- Sample holder for optical and electrical experiments
- Temperature Controller

#### **Options and Upgrades**

- 4.5 and 8 inch rotatable CF flanges available
- 4K Coldhead (0.1W @ 4.2K)
- 5.5K Coldhead (1W @ 10K)
- 450K High Temperature Interface
- 800K High Temperature Interface
- Turbo upgrade for faster cooldown times
- Custom temperature sensor configuration (please contact our sales staff
- Custom wiring configurations (please contact our sales staff)
- Window material upgrades (custom materials available)
- Sample holder upgrades (custom sample holders available)



The above picture shows a True UHV Closed Cycle Cryocooler



The above picture shows an instrumentation skirt with the electrical feedthroughs rotated 90 degrees upwards to allow for tight rotational



# **Cooling Technology**

DE-202	Closed Cycle Cryocooler
Refrigeration Type	Pneumatically Driven GM Cycle
Liquid Cryogen Usage	None, Cryogen Free
nperature*	

#### Tem

nperature*						
DE-202AI	< 10K - 350K					
DE-202PI	< 5.5K - 350K					
DE-202SI	< 4.2K - 350K					
With 800K Interface	(Base Temp + 2K) - 700K					
With 450K Interface	(Base Temp + 2K) - 450K					
Stability 0.1K						
*Based on bare cold head with a closed radiation shield, and no additional sources of experimental or parasitic heat load						

#### Sample Space

Diameter	Large Open Radiation Shield			
Height	Large Open Radiation Shield			
Sample Holder Attachment	1/4 - 28 screw			
Sample Holder	www.arscryo.com/Products/ SampleHolders.html			

#### **Chamber Interface**

Flanges	CF, ISO
Size	4.5", 6", 8", 10"

# Temperature Instrumentation and Control (Standard)

Heater	36 ohm Thermofoil Heater anchored to the coldtip
Control Sensor	Curve Matched Silicon Diode installed on the coldtip
Sample Sensor	Calibrated Silicon Diode with free length wires

Contact ARS for other options

#### **Instrumentation Access**

Instrumentation Skirt	True UHV Welded Stainless Steel
Instrumentation Ports	2
Instrumentation Wiring	Contact sales staff for options

# **Radiation Shield**

Material	Nickel Plated OFHC Copper
Attachment	Threaded
Optical Access	Open End Radiation shield terminates 0.125" short of cold tip (customer specified)


# **Cryostat Footprint**

Overall Length	466 mm (18.34 in)
Motor Housing Diameter	114 mm (4.5 in)
Rotational Clearance	200 mm (8 in) with "G" Configuration

Cryocooler Model		DE-202AB		DE-202A(T)B		DE-202PB		DE-202SB	
	Frequency	60 Hz	50 Hz	60 Hz	50 Hz	60 Hz	50 Hz	60 Hz	50 Hz
Base Temperature	•	<9K	<9K	<9K	<9K	<5.5K	<5.5K	<4.2K	<4.2K
Cooling Capacity*	4.2K	-	-	-	-	-	-	0.1W	0.08W
	10K	0.5W	0.4W	0.7W	0.56W	1W	0.8W	1.2W	1W
	20K	2.5W	2W	3.7W	3W	3.5W	2.8W	4W	3.2W
	77K	4W	3.2W	6W	4.8W	3.5W	2.8W	4W	3.2W
Radiation Shield C	ooling Capacity	10W	8W	15W	12W	10W	8W	10W	8W
Cooldown Time	20K	50 min	60 min	35 min	42 min	60 min	72 min	60 min	72 min
	Base Temperature	70 min	84 min	50 min	60 min	90 min	108 min	90 min	108 min
Compressor Model		ARS-	4HW	ARS-	4HW	ARS-	4HW	ARS-	4HW
Typical Maintenan	ce Cycle	12,000	hours	8,000	hours	12,000	) hours	12,000	hours

# True UHV — Low Power

# **CS202\*B Outline Drawing**

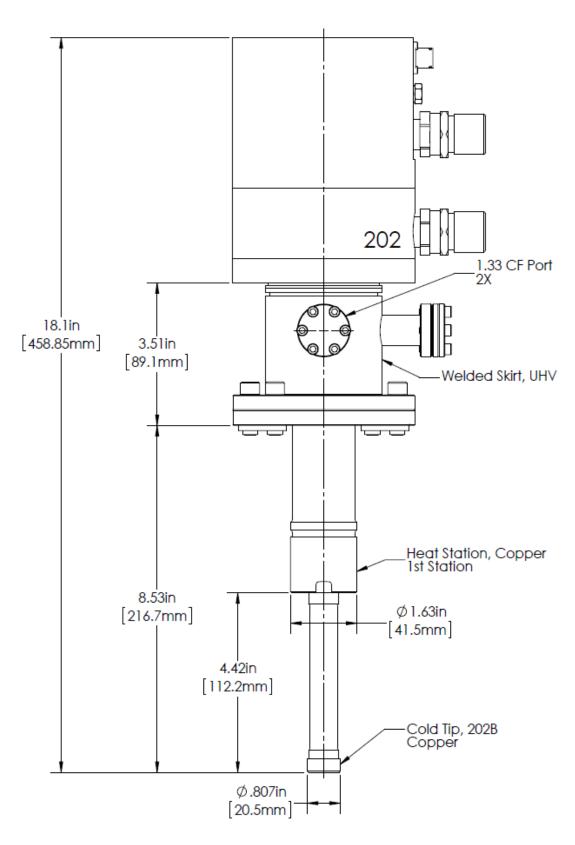




# True UHV — Low Power

# CoolPac



# **ARS-4HW Compressor**

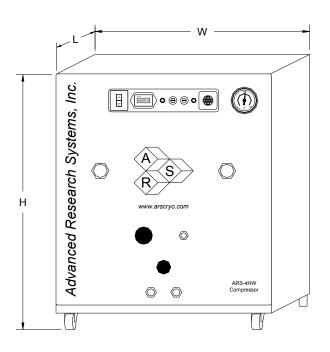

# **Compressor Model**

# ARS-4HW

	Frequency	60 Hz	50 Hz			
Standard Voltage	Min	208 V	190 V			
	Max	230 V	210 V			
Transformer Options	10%		220 V, 230 V			
	15%		240 V			
Power Usage	Power Usage Single Phase		3.0 kW			
Refrigerant Gas		99.999% Helium Gas, Pre-Charged				
Noise Level		60 dBA				
Ambient Temperature		12 - 40 C (54 - 104 F)				
Cooling Water	Consumption	2.3 L / min (0.6 Gal. / min)				
	Temperature	10 - 35 C (50–95 F)				
Connection		3/8 in. Swagelok Fitting				
Dimensions:	L	483 mm (19 in)				
	W	434 mm (17.1 in)				
	Н	516 mm (20.3 in)				
Weight		72 kg (160 lbs)				
Typical Maintenance Cyc	:le	12,000 hours				
	ion	CoolPac Comp				